	记录编号 FN: SMP-RD-A0002-R06	版本号 Ver: 2.1
	归档 Filing: filed immediately, permanently kept	生效日期 Ed: 2023.01.12

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多宁/DuoNing

动物细胞高性能培养基 High-Performance Culture Medium for Animal-Cells

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DN feed 4 & DN feed B4 are chemical defined (CD) highly efficient feed medium with no animal derived components, no protein, which are suitable for use in Fed-batch cultures using Chinese hamster ovary cells (CHO) in the development and production of therapeutic protein products. This product does not contain growth factors, hypoxanthine, thymidine or L-glutamine. It is suitable for the culture of different cell lines such as CHO-K1, CHO-S, DG44, CHOZN using GS system and DHFR screening system.

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DN feed 4 should be used together with DN feed B4, and added at the same time.

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The first replenishment is recommended in the middle to late stage of cell index growth, usually on day 3 when the cell density reaches $2.0\sim 3.0\times 10^6$ cells/mL or more, and the recommended feeding scheme is as follows, and the optimized ratio can be customized according to the cell line used:

补料时间 Feeding time	DN feed 4+DN feed B4
Day 3	3%+0.3%

	记录编号 FN: SMP-RD-A0002-R06	版本号 Ver: 2.1
	归档 Filing: filed immediately, permanently kept	生效日期 Ed: 2023.01.12

Day 5	3%+0.3%
Day 7	3%+0.3%
Day 9	3%+0.3%
Day 11	3%+0.3%

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DN feed 4 Liquid packaging: 2°C to 8°C, protect from light; validity period: 12 months.

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DN feed B4 Liquid packaging: 2°C to 8°C, protect from light; validity period: 12 months.

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